# P720 PERISTALTIC PUMPS

Instech's P720 peristaltic pumps deliver the accurate low flow rates of syringe pumps without the limits on delivered volume. They are ideal for animal IV infusion, tissue perfusion, and other low-flow laboratory applications.

The pump's analog circuitry has been carefully designed to minimize electromagnetic radiation; for this reason, the pump is often used for tissue perfusion even in the presence of sensitive intracellular recordings.

Instech offers this pump with four different motors; choose the motor and tube set size that best match your flow rate requirements. In general, use the standard P720 for flows between 0.05 and 1 mL/min, the lower flow P720/10K for rates between 5 and 100  $\mu$ L/min, the ultra low flow P720/37K for rates as low as 0.2 $\mu$ L/min, and then the higher-flow P720/66 for rates up to 18ml/min (1 L/hr). With a given tube set size, the pump performs best over a 10:1 flow control range.

Part No.	Description	Unit			
P720/37K	Ultra-low-flow peristaltic pump, tube sets not included	ea			
P720/10K	Very-low-flow peristaltic pump, tube sets not included	ea			
P720	Low-flow peristaltic pump, tube sets not included	ea			
P720/66	High-flow peristaltic pump, tube sets not included	ea			
P720/NCA	Fittings for straight-through tube sets	pair			
P720/RMC	Rod mounting clamp for P720 pumps	ea			
P720/KSK	Kapton® strip kit: 10 strips, 3 pcs grip paper, tool, clip	ea			
P720/BATT	9V lithium battery for P720 pump	ea			
S www.instechlabs.com/Pumps/peristaltic/					

S P E C I F I C A T I O N S

	P720/37K NEW	P720/10K	P720	P720/66		
Flow rates						
.015"ID tube .020"ID tube .031"ID tube .062"ID tube .093"ID tube	0.2 - 2.2 μL/min 0.4 - 4.6 0.9 - 8.7 3 - 30 6 - 58	0.8 - 7.5 μL/min 1.6 - 16 3 - 30 10 - 100 20 - 200	8 - 83 µL/min 18 - 180 35 - 350 125 - 1250 240 - 2400	0.07 - 0.7 mL/min 0.15 - 1.5 0.3 - 2.8 0.9 - 9 1.8 - 18		
Flow control range	~10:1	~10:1	~10:1	~10:1		
Accuracy	±5%	±5%	±5%	±5%		
Repeatability	±3%	±3%	±3%	±3%		
Motor gear ratio	36976:1	10683:1	900:1	66:1		
Gearhead type	spur	spur	spur	planetary		
Power source	Universal 15VDC 1A adapter, 2.5mm male plug, + tip Input: 100-240VAC, 50-60Hz; plugs for US, Europe, UK, Australia					
Backup battery	9V lithium	9V lithium	9V lithium	NA		
Backup battery life	~30 hours	~30 hours	~30 hours	NA		
Operating humidity	<80% (non-condensing)					
Dimensions	2.5 x 2.3 x 4.0" (6.4 x 5.7 x 10.2 cm)					
Weight	375g	375g	375g	375g		

For research and industrial applications only. Not approved or intended for human use.

### FEATURES

Analog speed control input ••••• for remote control or closedloop applications

Fine speed control

Coarse speed control



Rear view

P720/NCA

Fittings for straight-through tube sets are available when connectors are not desired

Snap-in tube sets with a wide range of tubing materials, sizes and connectors

Precision machined pump head with Swiss-made motor assembly for accuracy and repeatability

Kapton strip to improve accuracy ...... and tube set life

Battery backup for portability and protection during power failure (except P720/66 model)







# **P720 PERISTALTIC PUMPS**

#### Standard Tube Sets



A variety of tube sizes and connector types lets you tailor your P720 pump to your application. Tube sets typically last about one month under continuous operation.

New pumps do not include tube sets; please order them separately. The 12 standard configurations at right are stocked in moderate quantities for fast delivery. Tube sets are sold in packages of 5.

Part Number Inlet Tubing Outlet P720/TS-BS020SBS .020"ID silicone P720/TS-BS031SBS .031"ID silicone ŀn P720/TS-BS062SBS .062"ID silicone .062" barb .062" barb P720/TS-BS093SBS .093"ID silicone P720/TS-DBS031SDBS 2 x .031"ID silicone\* P720/TS-DBS062SDBS 2 x .062"ID silicone\* P720/TS-DBS031SDBSU 2 x .031"ID silicone, unbalanced1\* dual .062" barb dual .062" barb P720/TS-DBS062SDBSU 2 x .062"ID silicone, unbalanced1\* P720/TS-020S .020"ID silicone P720/TS-031S .031"ID silicone none<sup>2</sup> none<sup>2</sup> P720/TS-062S .062"ID silicone P720/TS-093S .093"ID silicone

<sup>1</sup>One tube intentionally shorter so that flow rate are lower. Use to maintain constant levels in tissue chambers. For details, see http://www.instechlabs.com/Pumps/peristaltic/levelcontrol.pl <sup>2</sup>Use with straight-through tube set fittings (P720/NCA).

(\$) www.instechlabs.com/Pumps/peristaltic/tubing.php

#### **Custom Tube Sets**

In addition to the standard silicone tube sets above, hundreds of custom co rations are possible. Create a part number as shown from the possible com tions and allow 4 to 5 days for the

e tu	ibe sets to be made to your spec	cifications.		þý—		<del>]-</del>			
	SIZE	015	020	031	062	093			
	Inner diameter	.015" (.38mm)	.020" (.51mm)	.031" (.79mm)	.062" (1.6mm)	.093" (2.4mm)			
	MATERIALS Use silicone with saline and most drugs. Use C-FLEX® for solutions containing fats, such as IV diets. Use VITON® for petroleum-based fluids.								
	S - Silicone	•	•	•	•	•			
	C - C-FLEX®		•	•	•				
	V - VITON®				•				
	CONNECTORS & CHANNELS								
	22 - 22ga tubing	•	•						
	20 - 20ga tubing	•	•	•					
	LL - male luer lock	•	•	•	•	•			
	FL - female luer lock	•	•	•	•	•			
	BS062" ID soft plastic barb	•	•	•	•	•			
	93093" ID soft plastic barb	•	•	•	•	•			
	D22 - dual channel 22ga tubing*	•	•						
	D20 - dual channel 20ga tubing*	•	•						
	DBS - dual channel .062" ID barb*		•	•	•				
	None (use with P720/NCA fittings)	•	•	•	•	•			
	C-FLEX is a reg trademark of Con	solidated Polymer Technologies. VI	ITON is a reg trademark of DuPor	nt Dow Elastomers. Tygon is a reg t	trademark of Saint Gobain Perforn	ance Plastics Corporation.			

\* DUAL TUBE SET WARNING: Dual channel tube sets, particularly the 062 size, place more stress on the pump motor than do single channel tube sets, which will shorten the life of the pump. Pumps with spur gearheads (P720/37K, P720/10K and P720) are more susceptible to this stress than are pumps with planetary gearheads (P720/66). In addition, dual 062 tube sets make contact with the sides of the pump rotor, which will cause those tube sets to wear faster than other sizes.

Instec	h_peri	staltic	pum	ps.qxp	111031



tube material

- outlet connector



LL Male luer loc Fits female luer

8



Fits .062" Tygon@



Fits PE50, .020" Tygon®

'n DBS Dual .062"ID barbs Fits .062" Tygon®

. 93 .093" ID barb



FL Female luer lock

Fits male luer lock



onfigu- Ibina- ons.		P	720/TS-F		
	020		031	062	
8mm)	.020" (.51	mm)	.031" (.79mm)	.062" (1.6mm)	

inlet connector-

tube ID